Using instructions for Adapters 1 x 180–250 ml

These using instructions are a supplement to the operating manual for the centrifuge and are not a substitute, so also read the operating manual before using the adapters for the first time. The current version of the operating manual and the safety instructions in your language can be found on the Internet at www.eppendorf.com.

1 Introduction

Adapters 1 x 180–250 ml are designed for swing-bucket rotor A-4-81 with rectangular buckets for Centrifuge 5810 / 5810 R. Bottles with a filling volume of 180–250 ml can be centrifuged in them. Conical tubes can also be centrifuged if an insertion form (ordered separately from the tube manufacturer) is used.


This adapter (gray) replaces the previous adapter (blue; order no. International 5825 720.008; order no. North America 022638921). You can use both adapters together, provided you follow the instructions below precisely.

2 Safety instructions and application limits


Risk of injury from asymmetrical loading of a rotor.

- ▶ Fill the rectangular buckets of the rotor with only two or four identical adapters for symmetrical loading.
- ▶ In doing this, be aware that the gray adapters differ significantly from the blue adapters in terms of weight.
- ▶ When using the gray and blue adapters together, only arrange them opposite one another: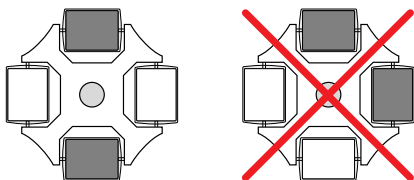

- ▶ Do not switch the module elements of the gray and the blue adapters, as this makes it impossible to tare the rotor:

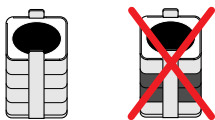Adapters with modules assembled correctly and incorrectly

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Eppendorf AG · Germany · Tel. +49 40 538 01-0 · E-Mail: eppendorf@eppendorf.com
Eppendorf North America, Inc. · USA · Toll free phone 800 645 3050 · E-Mail: info@eppendorf.com