

Instructions for use

Charger Shell for Charger Stand and Charger Carousel Series 4880

English (EN)

1 Exchanging the charger shell

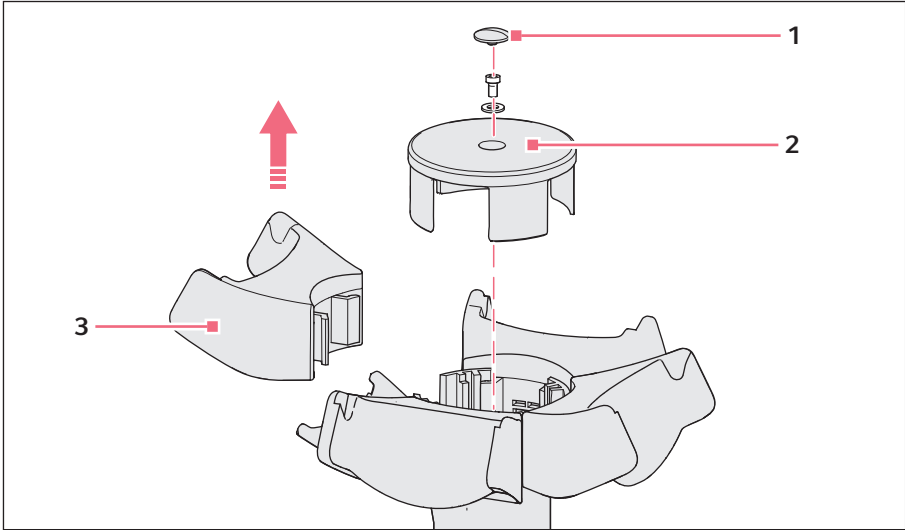


Fig. 1-1: Exchanging the charger shell

1 Sealing cap

3 Charger shell

2 Cover plate

To exchange a charger shell, proceed as follows:

1. Remove the charging plug of the power supply from the connector socket.
2. Remove the sealing cap (1).
3. Remove the screw on the cover plate (2). Use the 2.5 mm Allen key from the charger shell accessories for this purpose.
4. Store the screw and flat washer in a safe place for assembly.
5. Reach under the charger shells (3). Slightly tilt the charger shell on both sides and pull it upwards. Remove the charger shell together with the cover plate (2).

Instructions for use

Charger Shell for Charger Stand and Charger Carousel Series 4880
English (EN)

6. Insert new charger shells.

The sliding contacts of the charger shell must touch the gold-plated tracks of the charger carousel.

7. Assemble the charger carousel in reverse order.

8. Insert the charging plug of the power supply into the connector socket.

9. Check the charging function.

If you are connecting an Xplorer/Xplorer plus or a Multipette/Repeater (X)stream, *Device connected to power supply!* will appear on the display. When disconnecting the device, *Device disconnected from power supply!* will appear on the display. A flashing battery symbol is displayed during the charging process.

2 Ordering information

Order no. (International)	Order no. (North America)	Description
4880 000.000	4880000000	Charger stand 4880 for 1 Xplorer/Xplorer plus
4880 000.018	4880000018	for 1 Multipette/Repeater (X)stream
4880 000.026	4880000026	Charger carousel 4880 incl. power supply 4880 603.006 for 4 XplorerXplorer/Xplorer plus
4986 603.005	4986603005	Power supply with power plug adapters for Xplorer/Xplorer plus, Multipette/ Repeater (X) stream
4880 603.006	4880603006	for charger carousel series 4880
4880 601.003	4880601003	Charger shell for charger carousel series 4880 for Multipette/ Repeater (X)stream
4880 602.000	4880602000	for Xplorer/Xplorer plus

Your local distributor: www.eppendorf.com/contact

Eppendorf AG · 22331 Hamburg · Germany

eppendorf@eppendorf.com · www.eppendorf.com