eppendorf



New Brunswick™ Excella® E24/E24R Shaker

Operating manual

Copyright ©2018 Eppendorf AG, Germany. All rights reserved, including graphics and images. No part of this publication may be reproduced without the prior permission of the copyright owner.

Phillips® is a registered trademark of Phillips Screw Company in the United States and other countries.

Allen® is a registered trademark of Allen Manufacturing Company, USA.

Sticky Pad® is a registered trademark of American Covers, Inc., USA.

Eppendorf® and the Eppendorf Brand Design are registered trademarks of Eppendorf AG, Germany.

Excella® is a registered trademark of Eppendorf, Inc., USA.

New Brunswick™ is a trademark of Eppendorf AG, Germany.

Trademarks are not marked in all cases with TM or [®] in this manual.

Table of contents

1	Opera	ating instructions	. 5
	1.1	Using this manual	. 5
	1.2	Danger symbols and danger levels	. 5
		1.2.1 Danger symbols	. 5
		1.2.2 Danger levels	. 5
	1.3	Symbols used	. 6
	1.4	Abbreviations used	. 7
2	Safet	y	. 9
	2.1	Intended use	
	2.2	User profile	. 9
	2.3	Information on product liability	
	2.4	Application limits	
	2.5	Warnings for intended use	
		. J	
3	Produ	uct description	11
	3.1	Product overview	
	3.2	Inspection of boxes	
	3.3	Features	
	0.0	3.3.1 Operating modes	
		3.3.2 Visible/Audible user alarms	
		3.3.3 Data logging	
		3.3.4 Setpoint retention	
		3.3.5 Automatic restart	
		3.3.6 Drive interrupt	
		5.5.0 Bive interrupt	10
4	Instal	llation	15
•	4.1	Selecting the location	
	4.2	Unpacking the device	
	4.3	Utility requirements	
	4.4	Installation of platform	
	4.5	Flask clamp installation	
	4.6	Electrical connections	
	4.0	Liectrical connections	20
5	Oner	ation	21
3	5.1	Start/Stop	
	5.2	Continuous (untimed) run	
	5.3	Check setpoint	
	5.4	Timed functions	
	5.4 5.5		
	5.5	Alarm functions	
		5.5.1 Deactivating	
	г,	5.5.2 Reactivating	
	5.6	Temperature setpoint	
	5.7	Temperature offset calibration	
	5.8	Factory calibration	
	5.9	Power failure	
	5.10	Speed calibration	24

English (EN)

6	Opera 6.1	Control 6.1.1	panelStatus indicators	25 25
		6.1.2	LED display	
		6.1.3	Function indicators	
		6.1.4	User interface keys	26
7	Troub	leshooti	ng	27
	7.1		errors	
8	Maint	tenance .		29
	8.1		maintenance	
	8.2		q	
	8.3	Disinfed	ction/Decontamination	30
9	Techr	nical data	a	31
	9.1	Weight/	dimensions	31
		9.1.1	Device dimensions	31
		9.1.2	Platform dimensions	31
		9.1.3	Chamber dimensions	31
		9.1.4	Required space/footprint	31
	9.2	Applicat	tion parameters	32
		9.2.1	Speed	
		9.2.2	Capacity	
		9.2.3	Temperature	
	9.3	Power s	upply	32
10	Order	ing info	rmation	33
	10.1	Accesso	ries	33
11	Trans	port, sto	rage and disposal	35
	11.1	Disposa	I	35
	Index	· • • • • • • • • • • • • • • • • • • •		36
	Certif	icates		37

1 Operating instructions

1.1 Using this manual

- ▶ Read this operating manual completely before using the device for the first time. Also observe the instructions for use of the accessories.
- ▶ This operating manual is part of the product. Thus, it must always be easily accessible.
- ▶ Enclose this operating manual when transferring the device to third parties.
- ▶ You will find the current version of the operating manual for all available languages on our website www.eppendorf.com.

1.2 Danger symbols and danger levels

The safety instructions of this operating manual indicate the following danger symbols and danger levels:

1.2.1 Danger symbols

Hazard point	Heavy loads
Electric shock	Crushing
Biohazard	Burns
Explosion	Material damage

1.2.2 Danger levels

DANGER Will lead to severe injuries or death.	
WARNING May lead to severe injuries or death.	
CAUTION	May lead to light to moderate injuries.
NOTICE May lead to material damage.	

1.3 Symbols used

Example	Meaning
•	You are requested to perform an action.
1. 2.	Perform these actions in the sequence described.
•	List.
0	References useful information.

1.4 Abbreviations used

°C

Degree Celsius

cm

Centimeter

h

Hour

Hz

Hertz

kg

Kilogram

kWh

Kilowatt Hour

L

Liter

lb

Pound

in

Inch

m

Meter

min

Minute

mL

Milliliter

mm

Millimeter

Ы

Proportional/Integral

PPE

Personal Protective Equipment

rpm

Revolutions per Minute

Operating instructions New Brunswick™ Excella® E24/E24R Shaker English (EN)

s

Second

V

Volt

VA

Volt Ampere

VAC

Voltage in Alternating Current

2 Safety

2.1 Intended use

This device is exclusively intended for indoor use and for uniform movement and temperature control of biological solutions and cultures in reaction vessels.

2.2 User profile

The device may only be operated by trained lab personnel who have carefully read this operating manual and are familiar with the device functions.

2.3 Information on product liability

In the following cases, the designated protection of the device may be compromised.

The liability for the function of the device passes to the operator if:

- The device is not used in accordance with this operating manual.
- The device is used outside of the range of application described in the succeeding chapters.
- The device is used with accessories or consumables that were not approved by Eppendorf.
- Service or maintenance is completed on the device by people who are not authorized by Eppendorf.
- The owner has made unauthorized modifications to the device.

2.4 Application limits



DANGER! Explosion hazard

- ▶ Do not operate the device in areas where work is completed with explosive substances.
- ▶ Do not use this device to process any explosive or highly reactive substances.
- ▶ Do not use this device to process any substances which could create an explosive atmosphere.

Due to its design and the ambient conditions in its interior, the device is not suitable for use in potentially explosive atmospheres.

The device may only be used in a safe environment, e.g., the open atmosphere of a ventilated lab. The use of substances which may contribute to a potentially explosive atmosphere is not permitted. The final decision on risks associated with the use of such substances lies with the user.

2.5 Warnings for intended use



WARNING! Health risk due to poisonous, radioactive or aggressive chemicals

- ▶ Observe the national regulations for handling these substances as well as the material safety data sheets and manufacturer's application notes.
- ▶ Wear personal protective equipment (PPE).



CAUTION! Risk of injury and/or damage to unit

- ▶ A minimum load is required to safely operate the Excella 24/24R Shaker.
- ▶ Make sure the unit has a minimum load of 5.4 kg (12 lb) so that it can safely operate under the maximum speed (400 rpm). This includes the weight of the platform, flasks, and media.



CAUTION! Lack of safety due to incorrect accessories or spare parts

Accessories and spare parts that are not recommended by Eppendorf compromise the safety, function and precision of the device. Eppendorf cannot be held liable or accept any liability for damage resulting from the use of non-recommended accessories and spare parts.

▶ Only use accessories and original spare parts recommended by Eppendorf.



WARNING! Burns due to hot metal on the device and hot pistons

▶ Only touch the device and pistons when wearing protective gloves.



WARNING! Risk of crushing fingers with lid

- ▶ Do not reach between the lid and device, or into the lid locking mechanism, when opening and closing the device.
- ▶ Always fully open the lid so it cannot fall and close.



CAUTION! Risk to health due to lifting heavy loads

- ▶ Only lift the device with another person or using a suitable aid.
- ▶ Make sure to use a transport aid for transportation over long distances.



NOTICE! Damage due to overheating

- ▶ Do not place the device close to sources of heat (e.g., radiator, drying cabinet).
- ▶ Do not expose the device to direct sunlight.
- ▶ Ensure there is adequate distance to the wall and adjacent devices, on all sides of the device, in order to guarantee unobstructed air circulation.

3 Product description

3.1 Product overview

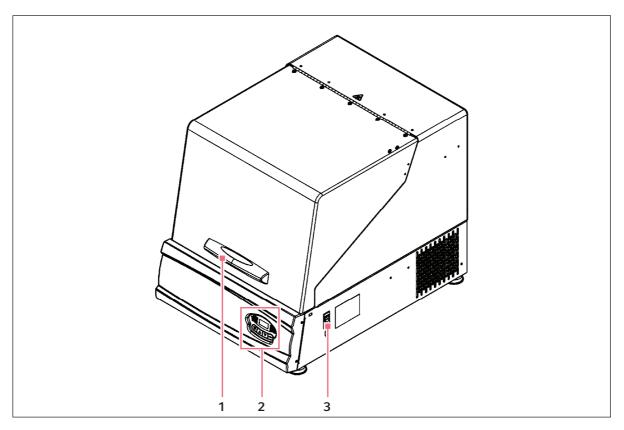


Fig. 3-1: Excella E24/24R shaker

1 Lid handle

2 User interface

3 ON/OFF mains/power switch

3.2 Inspection of boxes

After you have received your order, inspect the boxes carefully for any damage that may have occurred during shipping. Report any damage to the carrier and to your local Eppendorf distributor immediately.

3.3 Features

The New Brunswick E24/24R Shaker uses a UniCentric counter-balanced drive mechanism. It provide horizontal plane rotary motion in a 1.91 cm (¾ in) diameter circular orbit. A Proportional/Integral (PI) microprocessor controls the speed and temperature over the entire range.

The New Brunswick E24/24R Shaker operates from 7 °C above ambient to 60 °C. This range depends on relative humidity and other ambient factors. Ambient temperature is measured 1 m from the front of the unit.

Erlenmeyer flasks, 2.8 L Fernbach flasks, and a wide variety of tubes and plates can be accommodated using Eppendorf shaker accessories.

A wide variety of platforms can be used with the New Brunswick E24/24R Shaker:

- Universal platforms are the most flexible, providing hole patterns for flask clamps, test tube racks, and other accessories.
- Dedicated platforms are supplied with flask clamps attached; they are designed solely and expressly for this purpose.
- Test tube racks, microplate holders, and test tube rack holders are also available (a universal platform is needed for all test tube racks and holders).

For further information on these accessories (see Accessories on p. 33).

3.3.1 Operating modes

The New Brunswick E24/24R Shaker can be operated in the following modes:

- Continuously: at a set speed and temperature, until user intervention.
- **In a timed mode:** run at a set speed, time and temperature for a period of up to 99.9 h, after which the shaker automatically shuts off, while the temperature is maintained at its setpoint.

3.3.2 Visible/Audible user alarms

The New Brunswick E24/24R Shaker is equipped with visual and audible alarms for the following conditions:

- The end of a timed run
- Deviations from speed setpoint (5 min after lid is closed)
- Deviations from temperature setpoint (5 min after lid is closed)
- · Mains/power failure
- Lid open

3.3.3 Data logging

RS-232 connection provided for Data Logging.

3.3.4 Setpoint retention

All setpoints and operating status are retained in non-volatile memory.

3.3.5 Automatic restart

Automatic restart after power is restored, indicated by flashing display.

3.3.6 Drive interrupt

Drive interrupt shuts off power to the shaker when the lid opens. Acceleration/deceleration circuit prevents sudden starts and stops, minimizing both splashing and mechanical damage.

14

Product descriptionNew Brunswick™ Excella® E24/E24R Shaker English (EN)

4 Installation

4.1 Selecting the location



NOTICE! Damage due to overheating

- ▶ Do not place the device close to sources of heat (e.g., radiator, drying cabinet).
- ▶ Do not expose the device to direct sunlight.
- ▶ Ensure there is adequate distance to the wall and adjacent devices, on all sides of the device, in order to guarantee unobstructed air circulation.

Select the location according to the following criteria:

- Suitable mains/power connection according to the name plate
- · Stable, even, and resonance-free base
- Well-ventilated area and no direct sunlight to prevent additional temperature increases
- Ambient conditions of 10 °C 35 °C, 20 %RH 80 %RH non-condensing
- Able to accommodate 90.8 kg (200 lb)



Allow a minimum clearance of 7.6 cm (3 in) around the incubator shaker for ventilation, access to mains/power cord (rear panel), and access to mains/power switch and RS-232 port (right side).

4.2 Unpacking the device




Keep the packing material and transport securing device for later transport or storage.

- 1. Remove the packing material.
- 2. Remove the transport securing device.
- 3. Use the details included in the scope of delivery to check that the delivery is complete.
- 4. Check all parts for damage in transit. Contact Eppendorf Customer Service if parts are missing or transport damage is present.



Before placing the shaker on the bench, remove the 4 red plastic feet protectors.

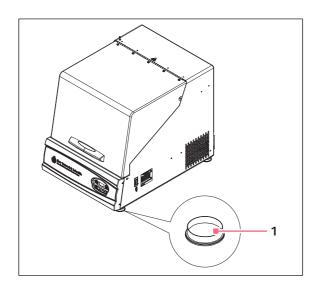

4.3 Utility requirements

The following utility requirements are needed for operation:

E24 mains/power supply

Utility	Requirement
Electricity	• 100 V, 50 Hz, 800 VA maximum
	• 100 V, 60 Hz, 800 VA maximum
	• 120 V, 60 Hz, 800 VA maximum
	• 230 V, 50 Hz, 800 VA maximum

E24R mains/power supply

Utility	Requirement
Electricity	 100 V, 50 Hz, 1500 VA maximum 100 V, 60 Hz, 1500 VA maximum 120 V, 60 Hz, 1500 VA maximum 230 V, 50 Hz, 1500 VA maximum

In all cases, voltage variations must not exceed ± 10 %.

4.4 Installation of platform



A platform with typical anticipated clamps and load must be installed prior to use.

The New Brunswick E24/24R Shaker is shipped with 4 Allen® head platform screws installed in the 4 bearing housing uprights. Before installing the platform, remove the screws and retain them for use.

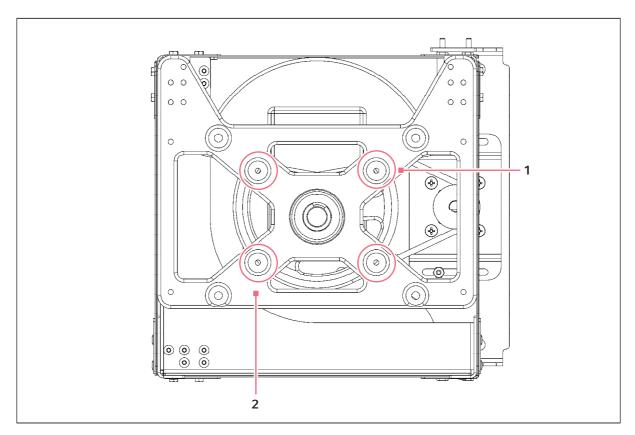


Fig. 4-1: Platform screw locations

1 Screws (4)

2 Sub-platform

- 1. Remove the 2 small plastic straps that secure the bearing housing in place for shipping.
- 2. Using the 7/32 in hex wrench provided, remove the 4 Allen head platform screws from the sub-platform. Save them for later use.
- 3. Place the selected platform on the sub-platform. Align the mounting holes of the platform with the platform screw locations in the sub-platform.
- 4. Insert the 4 Allen head platform screws previously removed. Tighten them with the 7/32 in hex wrench provided to secure the platform.

4.5 Flask clamp installation

Flask clamps purchased for use with the platform require installation. Clamps are installed by securing the clamp base to the platform with the correct type and number of screws. All clamps are shipped complete with hardware.



The New Brunswick E24/24R Shaker platform requires $10-24 \times 5/16$ in Phillips[®] head screws (which are supplied) to fasten flask clamps.

Clamps for 2.8 L and 2 L flasks are shipped with an extra girdle to keep the flasks in place. The girdle is an assembly of springs and sections of girdle tube. One girdle is already in place on the clamp, the other is packed separately.

To install these double girdle clamps:

- 1. Place the clamp on the platform, aligning its mounting holes with holes on the platform. Secure the clamp in place using the flask clamp screws provided (#S2116-3051, 10-24 x 5/16 in).
 - 3 different types of screws are shipped with the clamps, to identify the proper screws (see Fig. 4-2 on p. 18).
- 2. With the first girdle in place, as delivered, on the upper part of the clamp body, insert an empty flask into the clamp.
- 3. After making sure the sections of tubing are located between the clamp legs, roll the first girdle down the clamp legs as far as it can go. The tubing sections rest against the platform, and the springs are under the clamp base.
- 4. Place the second girdle around the upper portion of clamp body (just as the first girdle was initially). Make sure that its spring sections rest against the clamp legs, while its girdle tube sections sit against the flask, in between the clamp legs.





Fig. 4-2: Clamp fastener

Eppendorf flask clamps are used on various shaker platforms. Flat head screws of different lengths and thread pitch are used to secure the clamp. The following table identifies the proper screw for your shaker application by reference to the head style. Select the appropriate screws and set the others aside.

Description	Part number	Qty	Application
A	S2116-3051	25	5/16 in (7.9 mm) thick aluminum,
		pieces	phenolic and stainless steel platforms.
Flask clamp screws, $10-24 \times 5/16$ in,			



1 L and larger flask clamps are fastened with 5 screws.

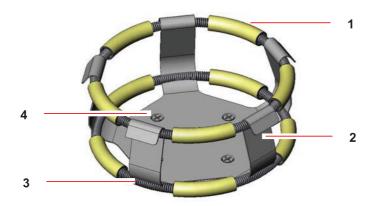


Fig. 4-3: Double girdle clamp installation

- 1 Upper girdle with girdle tubes
- 2 Clamp body (legs and base)
- 3 Lower girdle with girdle tubes
- 4 Clamp mounting holes (quantity: 5)



The upper girdle secures the flask within the clamp, and the bottom girdle keeps the flask from spinning.

4.6 Electrical connections

Please read the following warnings!



WARNING! Risk of electric shock and/or damage to unit

- Check that the voltage and frequency of your unit are compatible with mains/power supply.
- ▶ Set the mains/power switch to the OFF position.



WARNING! Risk of electric shock and/or damage to unit

▶ Use an earthed/grounded mains/power supply.



CAUTION! Risk of injury and/or damage to unit

- ▶ A minimum load is required to safely operate the Excella 24/24R Shaker.
- ▶ Make sure the unit has a minimum load of 5.4 kg (12 lb) so that it can safely operate under the maximum speed (400 rpm). This includes the weight of the platform, flasks, and media.
- ▶ Only after taking the preceding precautions, plug the mains/power cord into the mains/power supply.

5 Operation

5.1 Start/Stop



The New Brunswick E24/24R Shaker does not operate if the lid is open.

To initially start the shaker:

- 1. Close the lid.
- 2. Turn the **ON/OFF** switch to the **ON** position.

During start-up, the LED display indicates the model of your shaker.

3. To start or stop operation of the platform, press the **START/STOP** key on the keypad. When the shaker begins to operate, the LED display tracks the speed as it accelerates to the last setpoint entered.

5.2 Continuous (untimed) run

- 1. Press the **SELECT** key until the **RPM** indicator illuminates.
- 2. If the display indicates that the shaker is **OFF**, press the **START/STOP** key.
- 3. To enter the **SET** mode, press either the ▲ or ▼ key.

The **SET** indicator illuminates.

4. Set the speed by using the ▲ or ▼ key until the desired setpoint displays. Continued pressure the ▲ or ▼ key causes the setting to change more rapidly.



The setpoint can be changed during a run without stopping the shaker by following Steps 2 – 4 above. During speed changes, a visual alarm (flashing **RPM** indicator) flashes and an audible alarm sounds until the speed returns to within 5 rpm of the setpoint.

5.3 Check setpoint

- 1. Press the **SELECT** key until the desired indicator illuminates.
- 2. To enter the **SET** mode and display the current setpoint, briefly press either the ▲ or ▼key.



Holding the ▲ or ▼ key for more than 0.5 s causes the speed setpoint to change. Should this occur, resetting is necessary.

5.4 Timed functions

The shaker can be programmed to automatically stop after a preset time period of 0.1 h - 99.9 h. There must be power to the shaker in order to set the timer, although a timed run can be initiated while the shaker is either stopped or operating.

To set the timer:

- 1. Press the **SELECT** key until the **HRS** indicator illuminates.
- 2. To enter the **SET** mode and set the desired run time (between 0.1 h − 99.9 h, press either the ▲ or ▼ key.
 - 0

If the shaker is stopped, skip to Step 5. If the shaker is already running, continue to Step 3.

3. Press the **START/STOP** key.

The shaker stops and the display reads OFF.

4. Press the **START/STOP** key again.

The **TIME** indicator illuminates and the shaker starts the timed run.

If the shaker is stopped:

5. Press the **START/STOP** key.

The shaker starts the untimed mode.

6. Press the **START/STOP** key again.

The shaker stops and the display reads OFF.

7. Press the **START/STOP** key a third time.

The **TIME** indicator illuminates and the shaker starts the timed run.

To disable the visual alarm (flashing **TIME** indicator), press the **SELECT** key and change to any other function.

To cancel the timer without stopping the shaker, repeat Steps 1 and 2, and then immediately press the **START/STOP** key. The **TIME** indicator stops flashing and the display reads OFF.

5.5 Alarm functions

In addition to the visual alarm, the New Brunswick E24/24R Shaker has an audible alarm that is activated at predetermined times. It can be deactivated by using the **MUTE** function (see *Function indicators on p. 26*).

5.5.1 Deactivating

- 1. Press the **SELECT** key until the **MUTE** indicator illuminates.
- 2. To display **ON**, press the \triangle or ∇ key. Then press the **SELECT** key.

5.5.2 Reactivating

To reactivate the audible alarm:

- 1. Press the **SELECT** key until the **MUTE** indicator illuminates.
- 2. To display **OFF**, press the \triangle or ∇ key. Then press the **SELECT** key.

5.6 Temperature setpoint

Press the **SELECT** key until the function ${}^{\circ}$ C indicator illuminates. The temperature can be set from 5 ${}^{\circ}$ C above ambient temperature to 60 ${}^{\circ}$ C (non-refrigerated units) or from 4 ${}^{\circ}$ C – 60 ${}^{\circ}$ C (refrigerated units). Increasing or decreasing the setpoint is accomplished with the ${}^{\blacktriangle}$ or ${}^{\blacktriangledown}$ key.

During operation, if the temperature of the chamber is more than 1.0 °C higher or lower than the temperature setpoint, an alarm is triggered. This alarm consists of a flashing °C indicator and audible beep. The alarm automatically deactivates as the New Brunswick E24/24R Shaker achieves the set temperature.

5.7 Temperature offset calibration

The temperature probe and the temperature controller are calibrated together at the factory. The temperature probe measures the temperature of the air at the probe's location, near the heat exchanger return vent. The controller uses the probe input to adjust air temperature, up or down, to match the temperature setpoint.

Depending on various conditions within the chamber (e.g., flask placement and size, the heat produced by growing organisms, heat losses due to liquid evaporation from flasks), the display temperature may differ from temperatures within the flasks.

If you wish to have the temperature display ("Indicated Temperature") match the temperature at a given point, or match the average of a series of points within the chamber ("Actual Temperature"), proceed as follows:

- 1. Let the unit equilibrate at or near the desired temperature. Record the Indicated Temperature.
- 2. Record the Actual Temperature.
- 3. Calculate the temperature correction value: Actual Temperature Indicated Temperature = Temperature Correction Value.
- 4. Press the **SELECT** key until the °C indicator illuminates.
- Press the ▲ and ▼ keys simultaneously.
 The display indicates CAL.
- 6. Using the ▲ or ▼ key, enter the Temperature Correction Value calculated in Step 3.
- 7. To save the Temperature Correction Value to memory, press the ▲ and ▼ keys simultaneously.
 - 0

The °C indicator pulses rapidly to indicate it is not operating in the factory default mode.

5.8 Factory calibration

To return to the factory calibration:

- 1. Press the **SELECT** key until the °C indicator illuminates.
- 2. Simultaneously press the ▲ and ▼ keys. The display indicates CAL.
- 3. Using the ▲ or ▼ key, set the Temperature Correction Value to zero.
- 4. Simultaneously press the ▲ and ▼ keys. The rapid pulsing of the °C indicator stops.

5.9 Power failure

In the event of a mains/power failure, the New Brunswick E24/24R Shaker is equipped with an automatic restart function.

If the shaker was in operation prior to the power interruption, when power is restored the shaker begins to operate at its last entered setpoint. The LED display flashes and the audible alarm sounds, indicating that a mains/power failure has occurred. Press any key to stop the flashing in the display and the audible alarm.

5.10 Speed calibration

To calibrate the shaking speed:

- 1. Set the shaker to a speed that can easily be measured. If you are using a strobe, the minimum speed should be 250 rpm.
- 2. Compare the reading on the display to the measured reading.

If an adjustment is needed:

- 1. Press the **SELECT** key until the **RPM** indicator illuminates.
- 2. Press the ▲ and ▼ keys simultaneously. The display indicates CAL.
- 3. To change the displayed value to match the measured speed, press either the ▲ or ▼ key.
- 4. To save the adjustment, press the ▲ and ▼ keys simultaneously.
- 5. Using the mains/power switch, turn the shaker **OFF**, then turn it back **ON**.

6 Operating controls and function

6.1 Control panel

The control panel consists of the status indicators, LED display, function indicators, and the user interface keys.

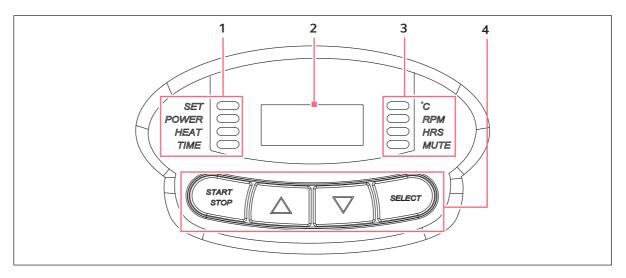


Fig. 6-1: Control panel

1 Status indicators

3 Function indicators

2 LED display

4 User interface keys

6.1.1 Status indicators

4 status indicator lights are located to the left of the LED display. They are:

SET Indicates that the shaker is in the **SET** mode, when setpoints are being

displayed and can be altered. This is activated by pressing the SELECT key or

by pressing the ▲ or ▼ key.

POWER Illuminates and blinks during power up or if power is interrupted during a run.

To turn off this indicator, press the **SELECT** key and change to another

function.

HEAT* Illuminates to indicate that the heater is on.

TIME Indicates that the timer is in operation. The shaker can be programmed to run

for a preset time from 0.1 h – 99.9 h. The timer can be disengaged without

stopping an ongoing run.

^{*}Not applicable for the E-2, E-5, and E-10

6.1.2 LED display

The digital display on the control panel is a three-digit LED display. During normal operation, the display indicates:

- Shaker status (On/Off)
- · Shaking speed
- · Chamber temperature*
- Setpoints
- · Hours remaining (in a timed run)
- Lid open ("LID")*

6.1.3 Function indicators

4 function indicator lights are located to the right of the LED display. They indicate the current parameter(s) being displayed:

°C* Interior chamber temperature. Can be set from 4 °C – 60 °C, when in **SET**

mode, using the ▲ or ▼ key. It indexes at 0.1 °C increments unless the key is

pressed for 4 s, after which it indexes more rapidly.

RPM Revolutions per minute. When in **SET** mode, use the ▲ or ▼ key to change the

speed. It indexes at 1 rpm increments unless the key is pressed for 4 s, after

which it indexes more rapidly.

HRS Time remaining in a timed run. Can be set from 0.1 h – 99.9 h, when in SET

mode, using the ▲ or ▼ key. It indexes at 0.1 h increments unless the key is

pressed for 4 s, after which it indexes more rapidly.

The countdown begins when the **START/STOP** key is pressed. If the **START/STOP** key is pressed, the shaking stops (but temperature is maintained) and

the timer pauses until the START/STOP key is pressed again.

When a timed run ends, the HRS indicator blinks. To turn off this indicator,

press the **SELECT** key and change to another function.

MUTE This feature is controlled by the **SELECT** key. When activated, the audible

alarm is muted, and remains so until is is reactivated. If MUTE is activated when the shaker is turned off using the ON/OFF switch, it remains engaged when the machine is powered up again. To activate (or deactivate) the MUTE function, press the SELECT key until the MUTE indicator illuminates. To turn the display ON or OFF, press the ▲ or ▼ key, and then press the SELECT key.

6.1.4 User interface keys

START/STOP Start or stop the shaker. It will also activate or stop the timer when a timed run

is desired.

SELECT Used to change the displayed parameter.

▲ or ▼ Used to adjust the setpoint of a displayed parameter up or down. They also

allow the user to enter the **SET** mode for setpoint changes.

^{*}Not applicable for the E2, E5, and E10

^{*}Not applicable for the E-2, E-5, and E-10

7 Troubleshooting7.1 General errors

Problem	Cause	Solution
Device does not run.	The power cable is loose.	Check the cable connections on the power connection and plug.
	The display is dark.	► Switch on the main power switch.
	The lid is open.	► Close the lid.
	The lid is closed.The open lid symbol is shown.	➤ The safety switch is defective. Contact your local Eppendorf service partner.
	The mains power switch is defective.	 Contact your local Eppendorf service partner.
Drive does not run at the set speed.	The device is overloaded.	Stop the drive.Reduce the sample weight.
	The device is loaded unevenly.	Stop the drive.Load the device uniformly.
Unusual operating noises	The device is loaded unevenly.	Stop the drive.Load the device uniformly.
	Foreign bodies or loose parts in the drive.	 Stop the drive and switch off the device. Contact your local Eppendorf service partner.
	Loose parts on the platform.	Stop the drive.Bolt loose parts on the platform.

8 Maintenance

8.1 Routine maintenance

No routine maintenance schedule is required for the New Brunswick E24/24R Shaker.

Clean the shaker occasionally using a cloth with conventional household (non-abrasive) cleaner.

To ensure proper air flow in and around the shaker, vacuum or sweep the area around the shaker to remove dust and other debris.

8.2 Cleaning



NOTICE! Risk of equipment damage

Use of abrasive or corrosive compounds can damage the shaker and void the warranty.

- ▶ Do not use abrasive or corrosive compounds to clean the shaker.
- 1. Routinely clean the exterior of the shaker by wiping it over with a soft cloth, moistened with soapy water
- 2. Rinse the soap from the cloth in clean water, and wipe the exterior surfaces again.

If there is biohazard contamination, see below.

8.3 Disinfection/Decontamination



WARNING! Risk of potential harm to personnel

▶ It is the responsibility of the user to carry out appropriate decontamination procedures if hazardous material is spilled on or inside the equipment. Before using any cleaning or decontamination method other than those suggested by the manufacturer, users should contact Eppendorf to ensure that the proposed method would not damage the equipment.



CAUTION! Risk of potential harm to personnel

- ▶ As a routine precaution, wear protective gloves.
- ▶ Be sure to adequately ventilate the work area as you disinfect, to avoid the formation of potentially explosive alcohol vapors.

Commercially available household bleach solutions, when diluted at a 1:10 ratio, are effective in routine decontamination of the shaker. The method for decontaminating a spill depends upon the nature of the spill.

- 1. Switch off the shaker. Unplug the shaker from the mains/power supply.
- 2. Spills involving fresh cultures or samples known to have low concentrations of biomass should be flooded with decontamination solution and soaked for 5 min before cleanup.
- 3. Spills involving samples with high concentrations of biomass, or involving organic matter, or occurring in areas warmer than room ambient temperature should be exposed to decontamination solution for at 1 h before cleanup.

9 Technical data

9.1 Weight/dimensions

9.1.1 Device dimensions

Width (external)	56 cm (22 in)
Height (external)	61 cm (24 in)
Height (with lid open)	101.9 cm (40 in)
Depth (external)	76 cm (30 in)
Weight (E24)	60 kg (133 lb)
Weight (E24R)	79 kg (175 lb)

9.1.2 Platform dimensions

Width	46 cm (18 in)
Depth	46 cm (18 in)
Weight	2.5 kg (5.5 lb)

9.1.3 Chamber dimensions

Width	51.7 cm (20 3/8 in)
Height	34.4 cm (13 9/16 in)
Depth	53.3 cm (21 in)

9.1.4 Required space/footprint

Width	68.6 cm (27 in)
Height	106.7 cm (42 in)
Depth	83.8 cm (33 in)

9.2 Application parameters

9.2.1 Speed

Speed (with 1.91 cm (3/4 in) orbit)	50 rpm – 400 rpm
Control accuracy	±1 rpm



Use of baffled flasks significantly reduces maximum speed for any shaker. We may be able to improve this maximum speed by using an alternative motor pulley. Contact Eppendorf Customer Service for more information.

9.2.2 Capacity

Maximum load	15.42 kg (34.00 lb)
Heater	Long-life, low-watt density resistance-type heater with high temperature thermostat

9.2.3 Temperature

Range (E24)	7 °C above ambient temperature to 60 °C	
Range (E24R)	15 °C below ambient temperature (minimum 4 °C) to 60 °C	
Accuracy	±0.1 °C at 37 °C	
Ambient operating environment	10 to 35 °C, 20 to 80 % relative humidity, non-condensing	
Altitude limit	2000 m	

9.3 Power supply

E24 mains/power supply

100 V, 50 Hz	800 VA
100 V, 60 Hz	800 VA
120 V, 60 Hz	800 VA
230 V, 50 Hz	800 VA

E24R mains/power supply

100 V, 50 Hz	1500 VA
100 V, 60 Hz	1500 VA
120 V, 60 Hz	1500 VA
230 V, 50 Hz	1500 VA

10 Ordering information

10.1 Accessories

Platforms	Part number
Universal platform	M1250-9902
125 mL Erlenmeyer flask dedicated platform*	M1194-9904
250 mL Erlenmeyer flask dedicated platform*	M1194-9905
500 mL Erlenmeyer flask dedicated platform*	M1194-9906
1 L Erlenmeyer flask dedicated platform*	M1194-9907
2 L Erlenmeyer flask dedicated platform*	M1194-9908
2.8 L Fernbach flask dedicated platform*	M1194-9932
Utility carrier with cushioned crossbars	M1194-9909
Utility tray with non-skid rubber surface	M1194-9910
Sticky Pad® platform	M1250-9903

^{*}dedicated platforms include flask clamps

Flask clamps	Part number
10 mL Erlenmeyer flask	ACE-10S
25 mL Erlenmeyer flask	M1190-9004
50 mL Erlenmeyer flask	M1190-9000
125 mL Erlenmeyer flask	M1190-9001
250 mL Erlenmeyer flask	M1190-9002
500 mL Erlenmeyer flask	M1190-9003
1 L Erlenmeyer flask	ACE-1000S
2 L Erlenmeyer flask	ACE-2000S
2.8 L Fernbach flask	ACSB-2800S

Racks and trays		Part number	Platform capacity
Adjustable angle test tube rack for	80 tube capacity	M1289-0110	4
tubes 8 mm – 11 mm diameter	60 tube capacity	M1289-0010	5
	48 tube capacity	M1289-0001	5
Adjustable angle test tube rack for	60 tube capacity	M1289-0200	4
tubes 12 mm – 15 mm diameter	44 tube capacity	M1289-0020	5
	34 tube capacity	M1289-0002	5
Adjustable angle test tube rack for	42 tube capacity	M1289-0300	4
tubes 15 mm – 18 mm diameter	31 tube capacity	M1289-0030	5
	24 tube capacity	M1289-0003	5
Adjustable angle test tube rack for	30 tube capacity	M1289-0400	4
tubes 18 mm – 21 mm diameter	23 tube capacity	M1289-0040	5
	18 tube capacity	M1289-0004	5
Adjustable angle test tube rack for tubes 22 mm – 26 mm diameter	22 tube capacity	M1289-0500	4
	16 tube capacity	M1289-0050	5
	13 tube capacity	M1289-0005	5
Adjustable angle test tube rack for tubes 26 mm – 30 mm diameter	20 tube capacity	M1289-0600	4
	16 tube capacity	M1289-0060	5
	12 tube capacity	M1289-0006	5
Microplate holder rack (stacked)	3 deep well or 9 standard	M1289-0700	8
Microplate holder rack (single layer)	5 deep well or standard	TTR-221	2
Angled test tube rack holder* for user-supplied test tube racks that are 10 mm – 13 mm (4 in – 5 in) wide and up to 38 mm (15 in) long.	NA	TTR-210	2
Angled test tube rack spacer bar* for use with TTR-210 to accommodate test tubes racks that are less than 13 mm (5 in) wide.	NA	TTR-215	NA

^{*}Universal platform required

11 Transport, storage and disposal

11.1 Disposal

In case the product is to be disposed of, the relevant legal regulations are to be observed.

Information on the disposal of electrical and electronic devices in the European Community:

Within the European Community, the disposal of electrical devices is regulated by national regulations based on EU Directive 2012/19/EU pertaining to waste electrical and electronic equipment (WEEE).

According to these regulations, any devices supplied after August 13, 2005, in the business-to-business sphere, to which this product is assigned, may no longer be disposed of in municipal or domestic waste. To document this, they have been marked with the following identification:



Because disposal regulations may differ from one country to another within the EU, please contact your supplier if necessary.

In Germany, this is mandatory from March 23, 2006. From this date, the manufacturer has to offer a suitable method of return for all devices supplied after August 13, 2005. For all devices supplied before August 13, 2005, the last user is responsible for the correct disposal.

Index	0
•	Operating speed32
A	
Accessories Flack clamps	P
Flask clamps	Platform 17
Platforms33	Power failure24
Racks and trays33	Power supply32
Actual temperature23	Product liability9
Alarm functions22	·
Audible alarm22	R
Auto restart function24	Requirements, utility
С	S
Control panel25	Speed calibration
D	Status indicators
Dimensions31	Symbols used6
Disposal35	т
_	Temperature offset calibration23
E 24 weight 21	Temperature setpoint23
E-24 weight	
E-24R weight31	U
F	User requirement9
Factory calibration24	Using this manual5
Flask clamp18	Utility requirements16
Function indicators26	
I	
Indicated temperature23	
Intended use9	
М	
Mains32	
Mains/power supply32	
Manual conventions6	
Mute alarm22	



Declaration of Conformity

The product named below fulfills the requirements of directives and standards listed. In the case of unauthorized modifications to the product or an unintended use this declaration becomes invalid.

Product name:

New Brunswick™ Excella® E24/E24R

including accessories

Product type:

Benchtop incubator shaker with optional refrigeration

Relevant directives / standards:

2014/35/EU: EN 61010-1, EN 61010-2-010, EN 61051-2-051

UL 61010-1, UL 61010A-2-10,

CAN/CSA C22.2 No. 61010-1, CAN/CSA C22.2 No. 1010.2.10

2014/30/EU: EN 61326-1, EN 55011

CISPR 11

EN 50419

Date: July 22, 2015

Management Board

Your local distributor: www.eppendorf.com/contact Eppendorf AG · 22331 Hamburg · Germany eppendorf@eppendorf.com

Eppendorf® and the Eppendorf logo are registered trademarks of Eppendorf AG, Germany. U.S. Design Patents are listed on www.eppendorf.com/ip. All rights reserved, incl. graphics and pictures. Copyright 2015 © by Eppendorf AG. ISO 9001 Certified

ISO 13485 Certified ISO 14001 Certified



Evaluate Your Manual

Give us your feedback. www.eppendorf.com/manualfeedback